


BUEHLER® EpoHeat™ EPOXY SYSTEM

- Heat Curing Epoxy With Excellent Clarity (Golden Color)
- Productive and Convenient: Reduce Time Spent in Tedious Weighing Operations by Making a Larger Batch Which Can be Poured for Up to 4 Hours Later
- Low Viscosity, Vacuum Impregnation and Heat Cure Results in Excellent Penetration of Pores, Cracks, and Other Fine Features
- Hard Epoxy With Low Shrinkage Gives Excellent Edge Retention
- Oven Cure Required



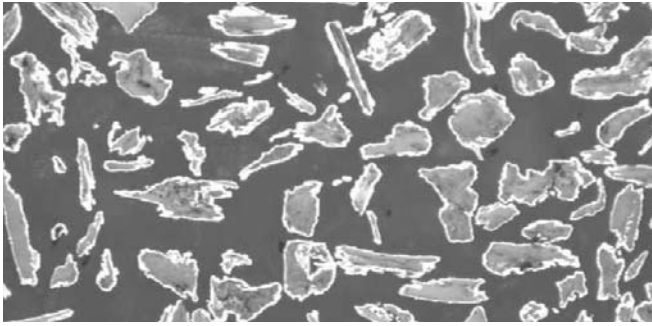

75Ni/C25 Nickel/Graphite powder mounted in EpoHeat 500X. Each individual flake was preserved during preparation and has excellent flatness and edge retention. Powder provided by Sulzer Metco (Canada) Inc.

- The luxury of a speedy process, before EpoHeat, was always at the sacrifice of another desired quality. Now achieve fast results with minimal shrinkage, excellent penetration and marvelous clarity with Buehler EpoHeat.
- Mixed viscosity approaches that of water during cure cycle, providing excellent penetration of pores, cracks, and other fine features
- Low viscosity allows EpoHeat to wet the substraight and fill voids that other epoxy systems cannot
- Heat curing assists in the impregnation of porous materials because trapped air molecules expand, rise to the surface and allow EpoHeat to penetrate the remaining void
- Extreme clarity means you can easily see into the mount for specimen reference and identification. Note color of mount is golden.
- Mix ratio by weight: 4 parts of Resin No. 20-8120 to 1 part of Hardener
- Can mix system twice a shift as mixed epoxy can sit for up to four hours without gelling
- Requires 90 minute oven cure at 131 °F (55 °C)
- May not fully cure, or cure correctly without oven cure
- Due to low viscosity for best results, use one piece molds
- Do not exceed oven temperatures of 167 °F (75 °C) as air bubbles and voids can develop if over exotherms
- Peak exotherm 295 °F (146 °C)
- 85 Shore D Hardness
- EpoHeat gives a 30% improvement in shrinkage over competitive systems resulting in improved preparation. Shrinkage Measurement Comparison: EpoHeat = 10µm, EpoThin = 12µm, Competitive System S = 15µm
- 1 year shelf life

EpoHeat Comparison to Other Buehler Castable Epoxy Systems

Product	Cure Time	Peak Temp	Shore D Hardness	HDT*	Recommended Mold
EpoKwick®	90 min	185 °F (85 °C)	82	130 °F (54 °C)	SamplKups®
EpoColor™	90 min	175 °F (79 °C)	82	110 °F (43 °C)	SamplKups®
EpoxiCure®	6 hours	82 °F (28 °C)	82	145 °F (63 °C)	Any
EpoHeat™	90 min	295 °F (146 °C)	85	170 °F (77 °C)	Reusable EPDM Mounting Cups
EpoThin®	9 hours	80 °F (27 °C)	78	110 °F (43 °C)	Any

*HDT (Heat Deflection Temperature) is a measure of distortion resistance under load at elevated temperature. HDT data shows EpoHeat will remain harder than our other epoxy systems during grinding/polishing.

Catalog Numbers

EpoHeat Resin 1qt. (0.95ℓ)20-8120-032
 EpoHeat Hardener 1 pt. (0.47ℓ)20-8122-016

Accessories

Release Agent with Swab** 2 oz. (0.03ℓ)20-8185-002
 Release Agent Refill ½ pt. (0.24ℓ)20-8185-008
 Release Agent Refill 1 pt. (0.47ℓ)20-8185-016
 Release Agent Refill 1 qt. (0.95ℓ)20-8185-032

** Used for mold release and other mounting applications

Recommended Molds

Resuable EPDM* Mounting Cups

Size	Quantity	Catalog Number
1" (2.5cm)	5	20-8181
1¼" (3.2cm)	5	20-8182
1½" (3.8cm)	5	20-8183
2" (5cm)	5	20-8184
2.2" x 1.2" x 0.9" (55mm x 30mm x 22mm)	1	20-8484
2.8" x 1.6" x 0.9" (70mm x 40mm x 22mm)	1	20-8485

* EPDM = Ethylene Propylene Diene Methylene Terpolymer

For a complete listing of Buehler Equipment and Consumables, please refer to our Buehler Equipment and Consumables Buyer's Guides. Buehler continuously makes product improvements; therefore, technical specifications are subject to change without notice.



BUEHLER
 BUEHLER LTD. – Worldwide Headquarters
 41 Waukegan Road • P.O. Box 1
 Lake Bluff, Illinois 60044-1699 USA
 Tel: 847/295-6500 • Fax: 847/295-7979
 Sales: 1/800/BUEHLER • 1/800/283-4537
 Web Site: <http://www.buehler.com>
 Email: info@buehler.com

BUEHLER GMBH - European and MESA Headquarters
 In der Steele 2 • 40599 Düsseldorf
 Postfach 16 03 55 • 40566 Düsseldorf
 Telefon: (49) 211 974100 • Telefax: (49) 211 9741079
 Web Site: <http://www.buehler-met.de>
 Email: info@buehler-met.de

BUEHLER FRANCE
 Téléphone: 0800 89 73 71
 Télécopie: 0800 88 05 27
 Web Site: <http://www.buehler.fr>
 Email: info@buehler.fr

BUEHLER UNITED KINGDOM
 Telephone: 0800 707 6273
 Fax: 0800 707 6274
 Web Site: <http://www.buehler.co.uk>
 Email: sales@buehler.co.uk

BUEHLER CANADA
 10 Carlow Court, Unit #2
 Whitby, Ontario L1N 9T7
 Telephone: (905) 430-4684 • Fax: (905) 430-4647
 Sales Telephone: 1-800-268-3593
 Email: info@buehler.ca

BUEHLER, ASIA-PACIFIC
 5/F Vogue Centre
 696 Castle Peak Road
 Lai Chi Kok, Kowloon
 Hong Kong, SAR, China
 Telephone: (852) 2307 0909
 Fax: (852) 2307 0233