



**Figure 1.** SEM view of Flat-Edge Filler (top, left). **Figure 2.** Example of a prepared mount using Flat-Edge Filler in EpoXide (bottom, left). **Figure 3.** Use of Flat-Edge Filler results in completely flat field from edge inward. Annealed 5140 alloy steel at 500x with Nital etch (right).

**BUEHLER®**

## Flat-Edge Filler

For Better Edge Retention with Castable Resins

Buehler proprietary Flat-Edge Filler is a mixture of fine globular ceramic oxides designed to improve castable resin abrasion resistance and edge retention.

The hardness of this filler is approximately 750 HV which is compatible with polishing of metals. This is much better than alumina filler, which at 1800 HV is too hard for grinding and polishing metals.

Buehler Flat-Edge Filler is chemically resistant and will not cause galvanic interactions during etching, as can happen when using certain plated metals for edge protection.

**Catalog No. 20-8196 0.45kg (1.0lbs)**

### Instructions for use

1. Mix up epoxy (Buehler EpoXide®, EpoKwik®, EpoColor®, or EpoThin®) as per instructions on epoxy container
2. Pour approximately ¼ of epoxy volume into a separate container and weigh (set aside the remaining ¾ volume of epoxy for use in filling up the mount later)
3. To the small ¼ volume of epoxy add 50% by weight, the flat-edge filler (one half part filler to one part epoxy)
4. Mix thoroughly together with stirring stick (approximately 1 minute)
5. Place the specimen into a mounting cup such as a SamplKup® and pour in filled epoxy mixture
6. Fill up the remainder of the mount with the normal unfilled epoxy
7. Allow to cure completely before preparing the sample as normal

**Note: Addition of the filler will slow down the cure cycle. If desired, place the mount in an oven at 80-100°F (27-38°C) to cure more quickly.**

*For additional information on our Flat-Edge Filler, please consult your nearest Buehler office. Buehler continuously makes product improvements; therefore, technical specifications are subject to change without notice.*

© 1999 BUEHLER LTD. Printed in U.S.A. .3M0799 FN01124



**BUEHLER**

**BUEHLER LTD. – Worldwide Headquarters**  
41 Waukegan Road • P.O. Box 1  
Lake Bluff, Illinois 60044-1699 USA  
Tel: 847/295-6500 • Fax: 847/295-7979  
Sales: 1/800/BUEHLER • 1/800/283-4537  
Web Site: <http://www.buehler.com>  
Email: [info@buehler.com](mailto:info@buehler.com)

**BUEHLER GMBH - European and MESA Headquarters**  
In der Steele 2 • 40599 Düsseldorf  
Postfach 16 03 55 • 40566 Düsseldorf  
Telefon: (49) 211 974100 • Telefax: (49) 211 9741079  
Web Site: <http://www.buehler-met.de>  
Email: [info@buehler-met.de](mailto:info@buehler-met.de)

**BUEHLER FRANCE**  
Téléphone: 0800 89 73 71  
Télécopie: 0800 88 05 27  
Web Site: <http://www.buehler.fr>  
Email: [info@buehler.fr](mailto:info@buehler.fr)

**BUEHLER UNITED KINGDOM**  
Telephone: 0800 707 6273  
Fax: 0800 707 6274  
Web Site: <http://www.buehler.co.uk>  
Email: [sales@buehler.co.uk](mailto:sales@buehler.co.uk)

**BUEHLER CANADA**  
10 Carlow Court, Unit #2  
Whitby, Ontario L1N 9T7  
Telephone: (905) 430-4684 • Fax: (905) 430-4647  
Sales Telephone: 1-800-268-3593  
Email: [info@buehler.ca](mailto:info@buehler.ca)

**BUEHLER, ASIA-PACIFIC**  
5/F Vogue Centre  
696 Castle Peak Road  
Lai Chi Kok, Kowloon  
Hong Kong, SAR, China  
Telephone: (852) 2307 0909  
Fax: (852) 2307 0233

**BUEHLER® Flat-Edge Filler**

